



SUMMARY


- **Duplex microcystin (MC) and cylindrospermopsin (CYN) direct toxin detection assay with quantitative results**
- **Field portable lysis method**
- **Can go from bloom to cell lysis and detection in 20 minutes**
- **Portable, low cost system**

SYSTEM FEATURES:

- **Multiplexed** – 6 or more (depending on assay)
- **Rapid** – 10 minutes assay plus 10 minutes lysis
- **Quantitative** – fluorescence immunoassay
- **Sensitive** – 0.5 ng/mL detection limit on MC
- **Robust** – field-ready technology
- **Versatile** – insensitive to debris in water sample


GEN 1 ASSAY FORMAT

Quantitative Duplex Immunoassay (MC and CYN)

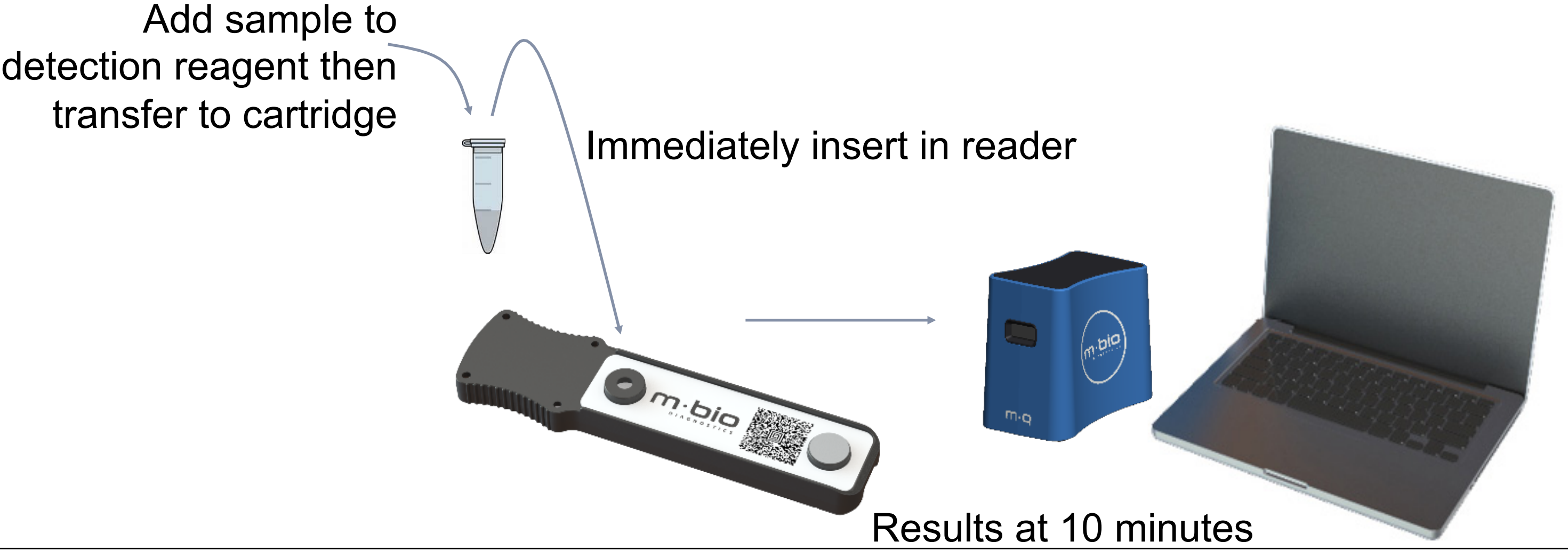


CALIBRATION / ANALYTICAL PERFORMANCE

- Factory calibrated which lasts for at least 6 months at room temperature and 12 months refrigerated
- Allows for the easy measurement of one or a few samples without the user measuring calibration curves
- Typical LOD is 0.5 µg/L for MC and 0.7 µg/L for CYN
- Typical standard curves for calibration are shown below

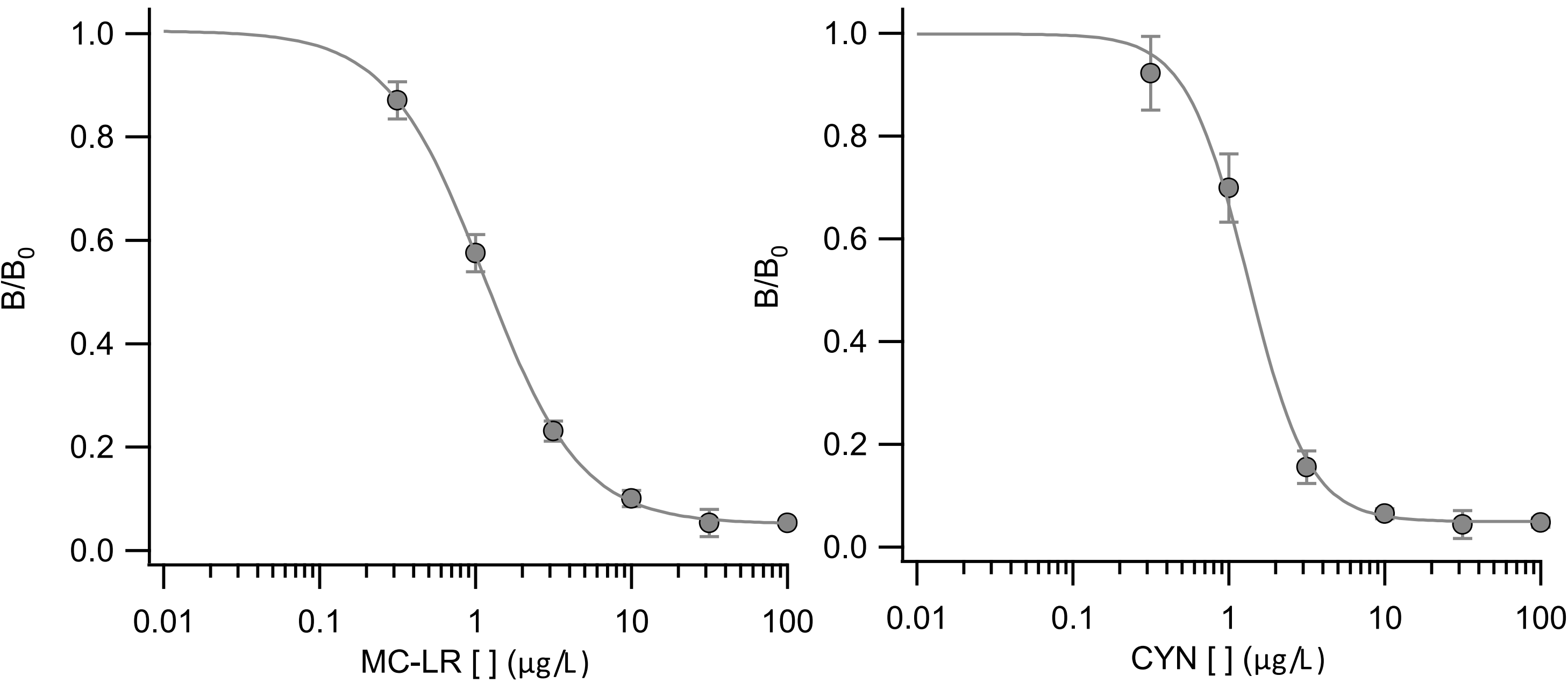

CONGENERS

- Goal is detection of total microcystin
- Proprietary monoclonal antibodies
 - Immunogen was an intact MC-LR conjugated to a carrier
 - Clone selected based on cross-reactivity against other congeners
- Dilution series performed with each of the congeners
- Certified Reference Materials
 - MC-LR, MC-RR, Nodularins, dM MC-LR
- Other Congeners
 - MC-LA, MC-LF, MC-YR, MC-LW

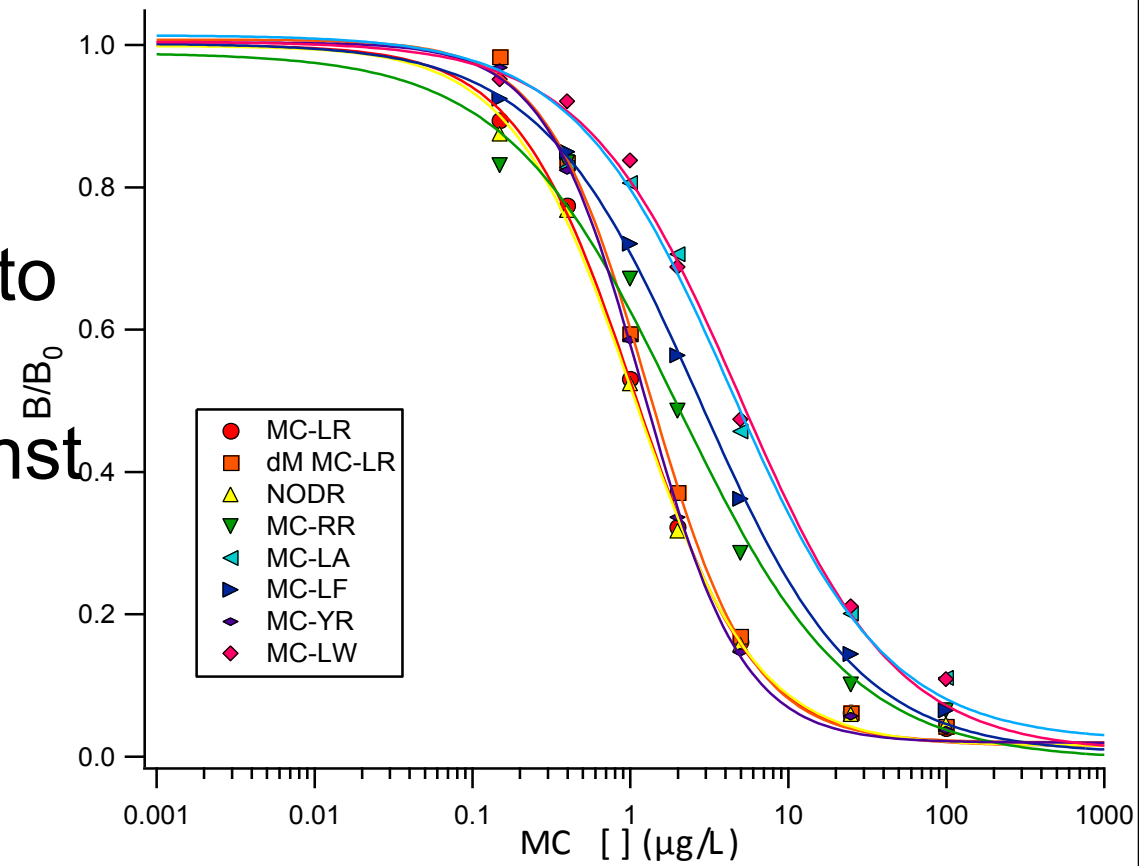

MQ Algae Lyse


- Rapid (10 minutes)
- Field portable, battery operated
- Effective (>95% lysis efficiency)
- Mechanical (bead beating) with a rotating whisk
- Measured toxin concentrations comparable to 3x freeze thaw cycles
- Cell lysis measured with microscopy

See Reference 1



MBio LIGHDECK[®] TECHNOLOGY

- Fluidic cartridge-based multiplexed immunoassays
- Novel planar waveguide-based illumination
- Fluorescence imaging detection
- Spatial multiplexing with microarray technology
- Quantitative output through proprietary software

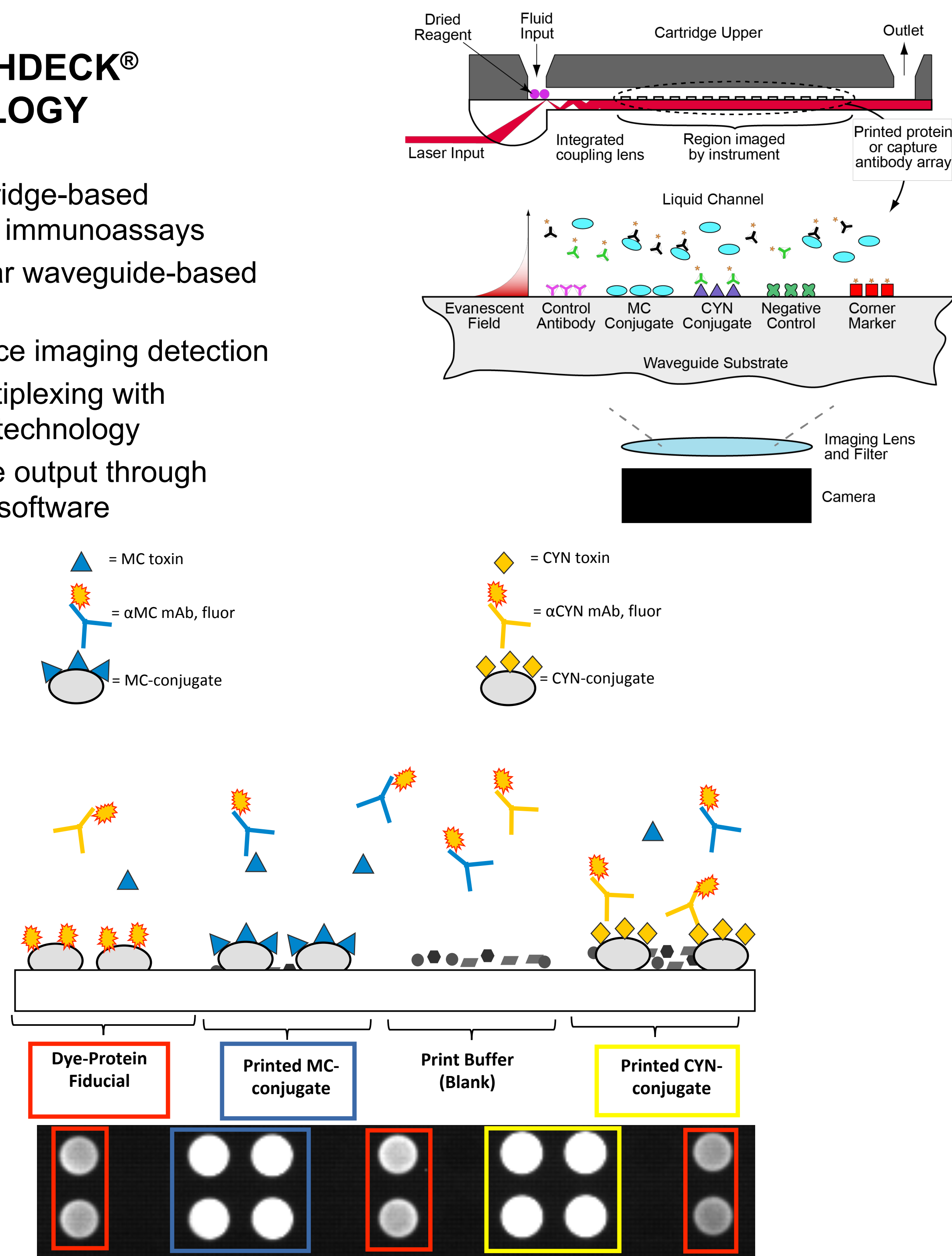

RESULTS ON NATURAL WATER SAMPLES

- 45 lake water samples were collected and tested on MBio platform and with a reference ADDA ELISA
 - 33 from Lake Erie
 - 12 from Colorado
- MBio results are the average of 3 measurements
- All samples were frozen and thawed 3x prior to measurement
- CYN was not detected by either ELISA or MBio in these samples
- Agreement between MBio and ADDA ELISA results for MC for Colorado and Western Basin of Lake Erie near Toledo
- In Sandusky Bay, Lake Erie, MBio reports lower concentrations of MC than ADDA ELISA, possibly due to congener cross reactivity
- Agreement of tested dilutions of CYN producing cell culture on MBio and ELISA

Colorado			
	Sample	MBio []	ADDA ELISA []
Active Bloom	Retention Pond A, Sample #1	16.54	20.25
	Retention Pond A, Sample #2	>80	>100
	Retention Pond B	<0.6	<0.15
No Bloom Visually Present	Aurora 1	<0.6	<0.15
	Aurora 2	<0.6	<0.15
	Aurora 3	<0.6	<0.15
	Boulder 1	<0.6	<0.15
	Boulder 2	<0.6	<0.15
	Westminster 1	<0.6	<0.15
	Westminster 2	<0.6	<0.15
	Westminster 3	<0.6	<0.15
	Westminster 4	<0.6	<0.15
		<0.6	<0.15

CYN Producing Cell Culture CS506		
Sample	MBio []	ELISA []
1:100 dilution	>2.7	2.10
1:200 dilution	1.50	1.67
1:300 dilution	0.93	1.12
1:400 dilution	0.63	0.76

Western Basin, Lake Erie near Toledo		
Sample	MBio []	ADDA ELISA []
GR1, Week 0	<0.6	0.19
GR1, Week 4	2.38	2.11
MB18 Week 4	6.42	6.66
MB20, Week 0	<0.6	0.22
7M, Week 0	>4.0	>5.0
7M, Week 1	<0.6	0.22
4P, Week 0	<0.6	<0.15
4P, Week 1	<0.6	0.35
CRIB, Week 1	<0.6	0.50
CRIB, Week 4	2.04	1.75
8M, Week 1	<0.6	0.29
8M, Week 4	4.52	6.38
Buoy (EW 5), Week 4	4.81	7.87

See Reference 1

Sandusky Bay, Lake Erie		
Sample	MBio []	ADDA ELISA []
Environment Canada Station 1163, Week 3	1.53	2.12
Environment Canada Station 1163, Week 5	<0.6	1.85
ODNR 1, Week 1	2.68	6.49
ODNR 1, Week 3	2.81	5.43
ODNR 2, Week 1	1.80	2.75
ODNR 2, Week 3	3.30	6.13
ODNR 2, Week 5	2.55	6.05
ODNR 4, Week 1	0.61	0.71
ODNR 4, Week 5	2.01	4.15
ODNR 6, Week 1	2.02	3.11
ODNR 6, Week 3	2.56	6.15
ODNR 6, Week 5	2.86	6.15
Sandusky Channel Bells, Week 1	1.03	1.25
Sandusky Channel Bells, Week 3	<0.6	0.27
Sandusky Channel Bells, Week 5	<0.6	0.49
Sandusky Buoy 2, Week 1	2.36	4.80
Sandusky Buoy 2, Week 3	2.15	5.24
Sandusky Buoy 2, Week 5	1.56	3.21
Edison Bridge, Week 3	2.56	6.91
Edison Bridge, Week 5	2.53	6.49

Comparison of lysis methods with cell cultures

Organism:	Microcystis aeruginosa		Cyl. ²
Identifier:	UTEX 2385	UTEX 2063	UTEX 942
Cell Conc. (cells/µL) x10 ⁴	2.1	1.6	1.2
Method	Percent Lysis		
3X Freeze-Thaw	62%	94%	76%
MBio Mechanical	84%	95%	99%

Comparison of lysis methods with natural water samples

Organism:	Aphanizomenon	Anabaena	Microcystis aeruginosa
Cell Conc. (cells/µL) x10 ⁴	1.3	0.3	24
Method	Percent Lysis		
3X Freeze-Thaw	99.8%	98.4%	98.7%
MBio Mechanical	99.2%	99.5%	99.9%

ACKNOWLEDGEMENTS

- We are grateful to our collaborators Greg Doucette, Tina Mikulski, Katrina Campbell, Christopher Elliott, Richard O'Kennedy, and Caroline Murphy.
- Samples were collected with assistance from Bowling Green State University, University of Toledo, Cities of Aurora, Boulder, Northglenn, Thornton, and Westminster, and Boulder Housing Partners.
- This material is based upon work supported by the NSF Grants No 1621951 and OCE-1440299. The views, opinions and/or findings expressed are those of the author and should not be interpreted as representing the official views or policies of the National Science Foundation or the U.S. Government.

REFERENCES

1. Bickman S, K. Campbell, C. Elliott, C. Murphy, R. O'Kennedy, P. Papst, M. Lochhead, An Innovative Portable Biosensor System for the Rapid Detection of Freshwater Cyanobacterial Algal Bloom Toxins, Environmental Science and Technology, 2018.
2. Reverte L., M Campas, BJ Yakes, J Deeds, P. Katikou, K. Kawatsu, M. Lochhead, C. Elliott, K. Campbell, Tetrodotoxin Detection in Puffer Fish by a Sensitive Planar Waveguide Immunosensor, Sensors and Actuators B: Chemical, 2017.
3. Devlin S, J. Meneely, B. Greer, C. Greef, M. Lochhead, C. Elliott, Next Generation Planar Waveguide Detection of Microcystins in Freshwater and Cyanobacterial extracts, utilizing a novel lysis method for portable sample preparation and analysis, Analytica Chimica Acta, 2013.
4. McNamee S, C. Elliott, B Greer, M. Lochhead, K. Campbell, Development of a Planar Waveguide Microarray for the Monitoring and Early Detection of Five Harmful Algal Toxins in Water and Cultures, Environmental Science and Technology, 2014.
5. Murphy C., E. Stack, S. Krivelo, D. McPartlin, B. Byrne, C. Greef, M. Lochhead, G. Husar, S. Devlin, C. Elliott, R. O'Kennedy, Detection of the cyanobacterial toxin, microcystin-LR, using a novel recombinant antibody-based optical-planar waveguide platform, Biosensors and Bioelectronics, 2014.